

**ROCK MAKER® Release Notes**

Last Updated: February 8, 2024

Version 3.18 Hotfix 3**February 8, 2024****Added:**

- A feature to configure the maximum age of retraining images sent to FORMULATRIX Scoring Server (FSS).
- Metadata for retraining images.

Fixed:

- An error when deleting an experiment with retraining images.
- Batchload import failure indicated by a pending status in the **Batchload Experiment Wizard**.
- Large log file creation when retraining images were disabled and the log level was verbose.
- Autoscoring malfunction when the ROCK MAKER Server detected corrupted images.

Version 3.18 Hotfix 1**June 23, 2023****Added:**

- The **Dispute** button in the **MARCO** panel in **Canvas** view.

Fixed:

- A failure to open an experiment in the ROCK MAKER Experiment Designer Edition.
- A failure to create new screens and open existing screens in the ROCK MAKER Experiment Designer Edition.
- Unresponsive or crashed **Explorer** in Windows 11.